R9000, RS9000 ***(HS-100D, HS-100D-Set)***

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Contents

1 Introduction	10
2 Safety Instructions	10
2.1 Description of Symbols	10
2.2 Instructions	10
3 Packing contents.....	11
4 Product description	12
4.1 Key features.....	12
4.2 General survey	12
4.3 Display description	12
4.4 Term description	13
5 Positioning the device and commissioning.....	13
6 Job description	13
6.1 Menu structure	13
6.2 Set speed	14
6.3 Set turn-off delay.....	14
6.4 Key lock	14
6.5 Storage function.....	15
7 Maintenance and cleaning.....	15
8 Technical data.....	15
9 Certificate.....	16

1 Introduction

Thank you for purchasing our **HS-D Overhead Stirrer**. You did a good choice buying a **Wisd** product.

Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you have to follow the safety instructions of the second chapter.

Please keep this operating manual in a save place for future reference and hand it out if you consign the device to a third person.

Witeg reserves to modify the capsule or make amendments for the improvement of quality and efficiency without prior notice.

2 Safety Instructions

2.1 Description of Symbols

**WARNING**

Information for avoiding injuries or fatal accidents.

**ATTENTION**

Safety instruction to avoid a damage of the product.

**ATTENTION**

Information about technical parameters.

**Environment**

Information about the operating environment.

**WARNING of explosions**

Safety instruction to avoid explosions.

2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Connect the device only to the suitable power supply. Please have a look to the identification plate.
- Use the device only in electric circuits with ground fault circuit interrupter.
- Check the power supply cord before you stick it into the electrical outlet.
- Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.

- Do not touch the device with wet hands during the usage because this might cause an electrical shock.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.

- Use this device only if you read the whole safety instructions.
- This device should only be used by persons who are familiarized with safety instructions made for laboratories.
- Pay attention that no liquid reaches the controller or inward of the capsule.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective work clothing during the usage.
- Pull out the power cable if you didn't want to use the device for a longer time.

- Do not use this device in areas with potentially explosive atmosphere.

- Make sure that the device can not slip away.
- The device could get damaged if the vibrations are too strong.
- To avoid heat congestions or fires make sure that there is enough space for air circulations.
- Avoid direct solar radiation.
- Do not use the device near highly combustible materials.
- Do not set the device outdoors.
- It is superiorly to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- To set the device in wet or dusty places could cause overheating, short circuits or fires.

- All circuits and electrotechnical components are **witeg** owned technology. In fact of this it is necessary that only **witeg** authorized personal repair the device. Contraventions exclude the warranty.

3 Packing contents

quantity	DH.WOS01015	DH.WOS01025	DH.WOS01016	DH.WOS01026
1	HS-30D	HS-100D	HS-30D-Set	HS-100D-Set
1	Allen key			
1	body holder			
1	chuck key			
1	operating manual			
1			stand rod ST110	stand rod ST120
1			clamp CL200	
1			impeller PL010	impeller PL015

4 Product description

4.1 Key features

The **HS-D Overhead Stirrer** is designed for chemical, biological, medical and pharmaceutical applications.


Key features are as follows:

- Jog-Shuttle knob for easy and fast handling.
- Up to 3000 rpm. Max. viscosity: 50.000 mpas.
- Constant speed in spite of viscosity fluctuation.
- Motor provided with overload and overheat protection.
- Stepless variable adjustability of every stirrer concerning to the continuous bored chuck.
- Power bar graph and speed showed on illuminated LCD.
- Smooth increase of speed for higher safety.
- Storage functions for speed value.

4.2 General survey

4.3 Display description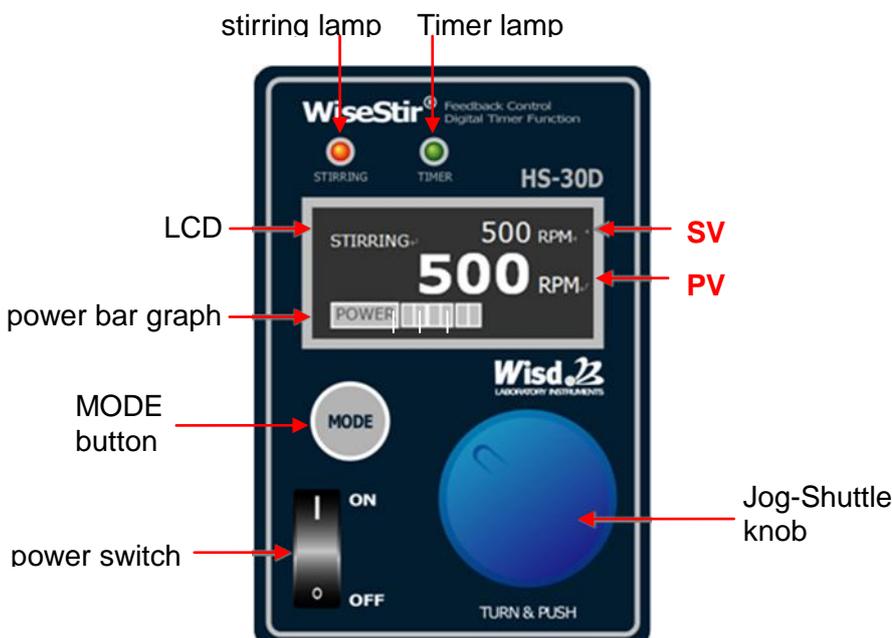

4.4 Term description

SV: set value, upper line in display, in key lock only PVs

PV: present value, lower line in display, in key lock upper line too

MODE button

- For switching between the different menu items.

Jog-Shuttle knob

- Change SVs by turning.
- Push the knob for start and stop or probate and store set values.

Main switch

- On / off switch.

Timer lamp

- Is blinking in menu items for timer modulation.
- Shines permanent in key lock if delay time is running.

Stirring lamp

- Is blinking in menu items for speed modulation.
- Shines permanent during the stirring process.

5 Positioning the device and commissioning


1. Follow especially the safety instructions marked with following symbol: 
2. Remove the package and perhaps soiling carefully.
3. If there are any unexpected transport damages, please inform your forwarder and your customer service.
4. Attach the clamp to the tripod rod.
5. Stick the device at the clamp and tighten the screw.
6. Check that each connection fits close.
7. Fix the stirrer in the chuck with included chuck key.
8. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatised.
9. Make sure that you follow all safety instructions.
10. Connect the device with the power outlet.
11. Switch the device on and start the process.

6 Job description

6.1 Menu structure

The **HS-D Overhead Stirrer** proffers you the genial combination of MODE button and Jog-Shuttle knob. That allows a simple and comfortable usage.

The menus structure is well arranged and pictured as follows:

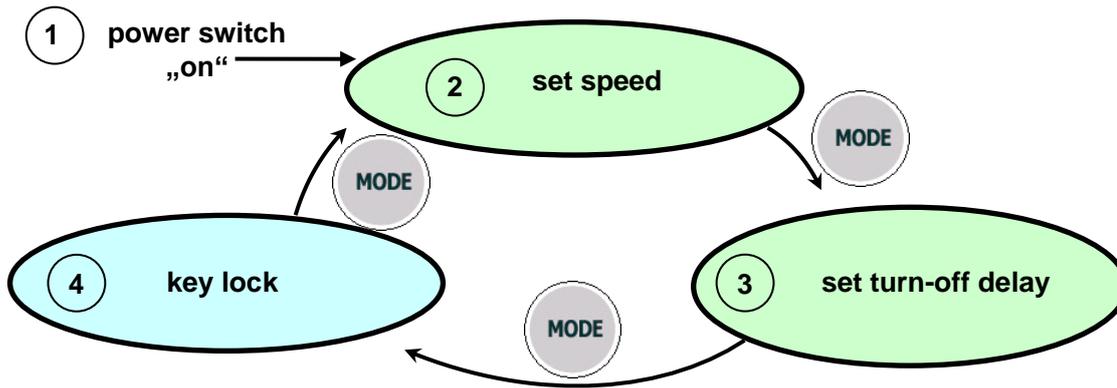6.2 Set speed

- First menu item after switching the device on is the menu “set speed”. The stirring lamp is blinking. You can recognize the menu at „RPM“ in SV- und PV-line.
- If there is showed “O RPM” in the SV line, there is no value stored and you can start the stirring process by turning the Jog-Shuttle knob clockwise.
- If “STIRRING” is blinking on the LCD there is a speed value stored. The saved value is showed in the SV-line and you can start the stirring process by pushing the Jog-Shuttle knob.
- In both cases is the speed during the stirring process variable. To change the SV you only have to turn the Jog-Shuttle knob clockwise to increase the speed and counter-clockwise to reduce it.
- To stop the stirring process manually, you have to push the Jog-Shuttle knob again or decrease the speed till “0” by turning the Jog-Shuttle knob counter-clockwise.
- The set speed is increasing gradually, so it could takes some time till the SV is climaxed.

6.3 Set turn-off delay

- The delay timer serves as automatically stop function. It turns off the process after elapsing the timeframe you set.
- To reach the timer menu you have to push the MODE button in the menu item “set speed”. The timer lamp is blinking.
- By renewed using of the MODE button you reach the key lock without delay time setting.
- The delay time is adjustable from one minute till over 99 hours (00:01 - 99:59) in one minute steps.
- Turn the Jog-Shuttle knob to set your delay time. Push the Jog-Shuttle if you reached the desired value to confirm it. With the confirmation the timeframe is elapsing
- If you want to change the value after confirmation, you have to push the Jog shuttle once again. The PV initiates blinking and you can modify the value again.
- To confirm the value push the Jog-Shuttle knob yet again.

6.4 Key lock

- To avoid accidental changing of the set values, push the MODE button in the menu item “set delay time”.
- In this menu item can you see the actually speed in the SV-line and the remaining time of delay function in the PV-line. The function of the Jog-Shuttle knob is deactivated.
- Press the MODE button once more to leave the key lock. Than you are in the menu item “set speed”.

6.5 Storage function

The **HS-D Overhead Stirrer** includes a storage function for speed and delay values. That is practical if you want to work with the same parameters often. Then you need not to set them in every usage.

- For storing values you only need to hold the Jog-Shuttle knob pushed for two seconds. Successful storing is displayed by blinking of the SV-value for three times.
- If you saved a speed value, "STIRRING" is blinking after the device's restart.
- If every stored parameter is erasable. Turn the value with the Jug-Shuttle till "0" and then hold the knob pressed for two seconds.
- If you want to store a new value it is not necessary to erase the old one.

7 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device.
- Pull the power plug out of the electrical outlet.
- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- Only use original spare parts.

8 Technical data

		HS-30D	HS-100D
Stirring Capacity (H₂O)		max. 10 Lit.	max.20 Lit.
Max. Viscosity (mPas)		10.000	50.000
Motor Rating (Input / Output)		83 / 75 W	130 / 120 W
Maximum Torque (Ncm)		85	145
Max. Shaft Diameter (mm)		8	
Speed Range (rpm)		200 – 3.000	
Controller		Digital Feedback Controller with Jog Shuttle knob (Turn + Push)	
Display		Digital LCD Display (RPM and Power bar-graph)	
Operating Environment	Ambient Temp. Range (°C)	0 ~ 40	
	Max. Permissible Humidity (%)	80	
Safety Device		Overload & Heat Protector for motor	
Dimension (wxdxh mm)		90x145x180	
Weight (kg)		4,5	5,3
Others		Smooth Start and Change of the Stirring Speed Storage Function	
Power Supply		AC 230 V, 1N~, 50/60 Hz	

9 Certificate

Certificate

for

The quality and all features were checked by the manufacturer before the shipment.

We grant from date of purchase

two years guaranty.

This certificate excludes damages by natural disasters or incorrect usages by the costumer.

Please look on your account and complete following table:

Article	HS-D Overhead Stirrer
Typ	R9000, RS9000 (HS-100D, HS-100D-Set)
Serialno.	
Date	

witeg Labortechnik GmbH

Am Bildacker 16
D-97877 Wertheim
TEL: +49-9342-9301-0
FAX: +49-9342-9301-77
Email: info@witeg.de
www.witeg.de